

HumanKine[®] IL-23 100x More Potent For Induction of Human Th17 Cells

INTRODUCTION

Cytokines are a group of proteins and polypeptides that organisms use as signaling molecules. Most cytokines are glycoproteins less than 30 kDa in size and bind to specific, high-affinity cell surface receptors. Due to their central role in the immune system, cytokines are involved in a variety of immunological, inflammatory and infectious diseases and widely used in research, diagnostics and therapeutics. Currently, these proteins are predominantly produced in non-human cells (e.g. E. coli, SF9, CHO) and therefore lack authenticity due to the absence of physiologically relevant glycosylation. In addition, a number of important cytokines are not commercially available due to inadequate proteolytic processing, protein folding or other post-translational modifications that do not occur in the non-human cell expression systems. Proteintech has developed an efficient human-cell based technology, HumaXpress[®], for the scalable production of human cytokines. The company is expanding the range of tag-free produced cytokines, including difficult-to-express members of the TGF β superfamily. HumanKine cytokines are preferred reagents for stem cell, cancer, inflammation research, and antibody development.

IL-23

Currently, commercially available recombinant IL-23 cytokine is produced as a heterodimeric or fusion protein from an insect cell expression system. Proteintech has produced HumanKine IL-23 in a stable cell culture of engineered human HEK293 cells. The protein is expressed as a disulfide-linked dimer of 55 kD and, due to the scalability of the stable culture, can be cost-effectively produced (Fig. 1).

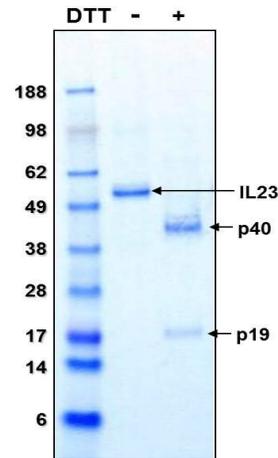


Fig. 1 Highly purified recombinant human Humankine IL-23 from human cells which is glycosylated and dimeric

IL-17 producing CD4⁺ T cells (Th17 cells) have been identified as a unique subset of T helper cells that develop along a pathway that is distinct from the Th1 and Th2-cell differentiation pathways. This finding has provided exciting new insights into immunoregulation, host defense and the pathogenesis of autoimmune diseases. Recently it has been shown that TGF β 1, IL-1 β , IL-6, and IL-23 are important in driving human Th17 differentiation. The bioactivities of IL-23 from human and insect cells were first determined by the dose-dependent secretion of IL-17 from mouse splenocytes activated with 10 ng/ml PMA, which shows that HumanKine IL-23 is tenfold more active (Fig. 2). The activities were further assayed with human CD4⁺ cells which were isolated from a healthy donor and stimulated with 10 μ g/ml plate bound anti-CD3 and 10 μ g/ml soluble anti-CD28 in the presence of Th17 polarizing cytokines. After 5 days supernatants were harvested for measurement of IL-17 by ELISA. The results show that HumanKine IL-23 is 100-fold more potent for inducing IL-17 secretion

in two independent studies; maximum induction was achieved with 0.1 ng/ml HumanKine IL-23 vs. 10 ng/ml with insect cell-produced IL-23 (Fig. 3). These results demonstrate that authentic human cell expressed cytokines can induce Th17 cell differentiation at physiologically relevant concentrations and may lead to a more accurate scientific understanding of the human biological process.

A rapidly expanding range of HumanKine cytokines are available from Proteintech Inc. The proteins are manufactured to high quality standards and provide high biological activity, lot-lot consistency and low endotoxin levels. The specific products discussed here, HumanKine IL-23, HumanKine TGF β 1, HumanKine IL-1 β , and HumanKine IL-6 are available for purchase from Proteintech.

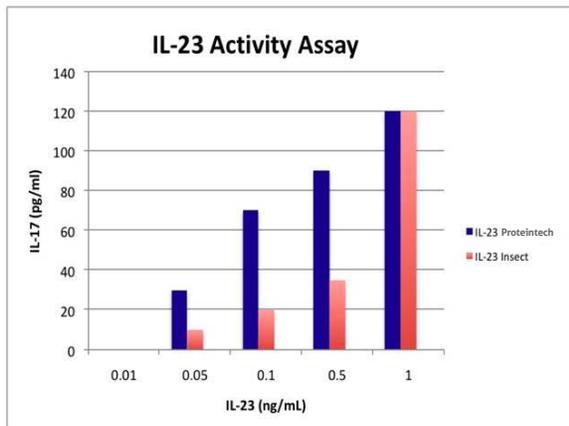


Fig. 2 IL-23 activities determined by the dose-dependent secretion of IL-17 from mouse splenocytes activated with 10 ng/ml PMA.

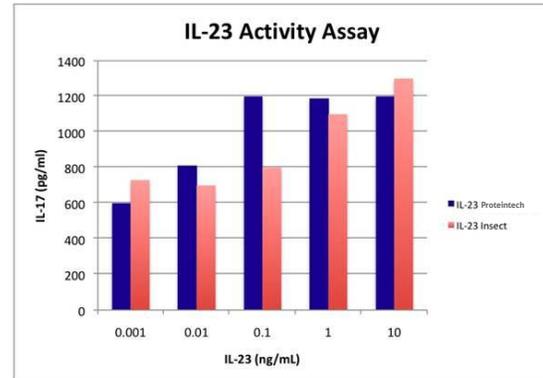


Fig. 3 IL-23 activities determined by the dose-dependent secretion of IL-17 from human CD4+T cells stimulated with 2 mg/ml plate bound anti-CD3 and 1 mg/ml soluble anti-CD28 in the presence of Th17 polarizing cytokines.